



Combining the Power of Mass Spectrometry with Ion Chromatography through a Single Software Solution

*Environmental Monitoring Conference, Crystal City, VA USA
August 1st – 5th, 2022*

Presented by

Jay Gandhi, *PhD*

Disclaimer of Endorsement

Mention of or referral to commercial products or services, and/or links to non-EPA sites does not imply official EPA endorsement of or responsibility for the opinions, ideas, data, or products presented at those locations, or guarantee the validity of the information provided. Mention of commercial products/services on non-EPA websites is provided solely as a pointer to information on topics related to environmental protection that may be useful to EPA staff and the public.

Today we will talk about.....

- **Perchlorate**
 - Background and History
 - Ion Chromatography methods
 - Ion Chromatography with Hyphenated Methods
- **What else can we analyze along with Perchlorate?**
- **Summary**


Perchlorate Background

- Used as rocket propellant
- Used in electroplating industry
- Used in fireworks
- There are also evidence of "*naturally occurring Perchlorate – Chilean fertilizer*"

Why Analyze?


- It is persistent in environment
- It is believed to inhibit Iodine uptake in thyroid gland causing *hypo* thyroidism

Perchlorate Background

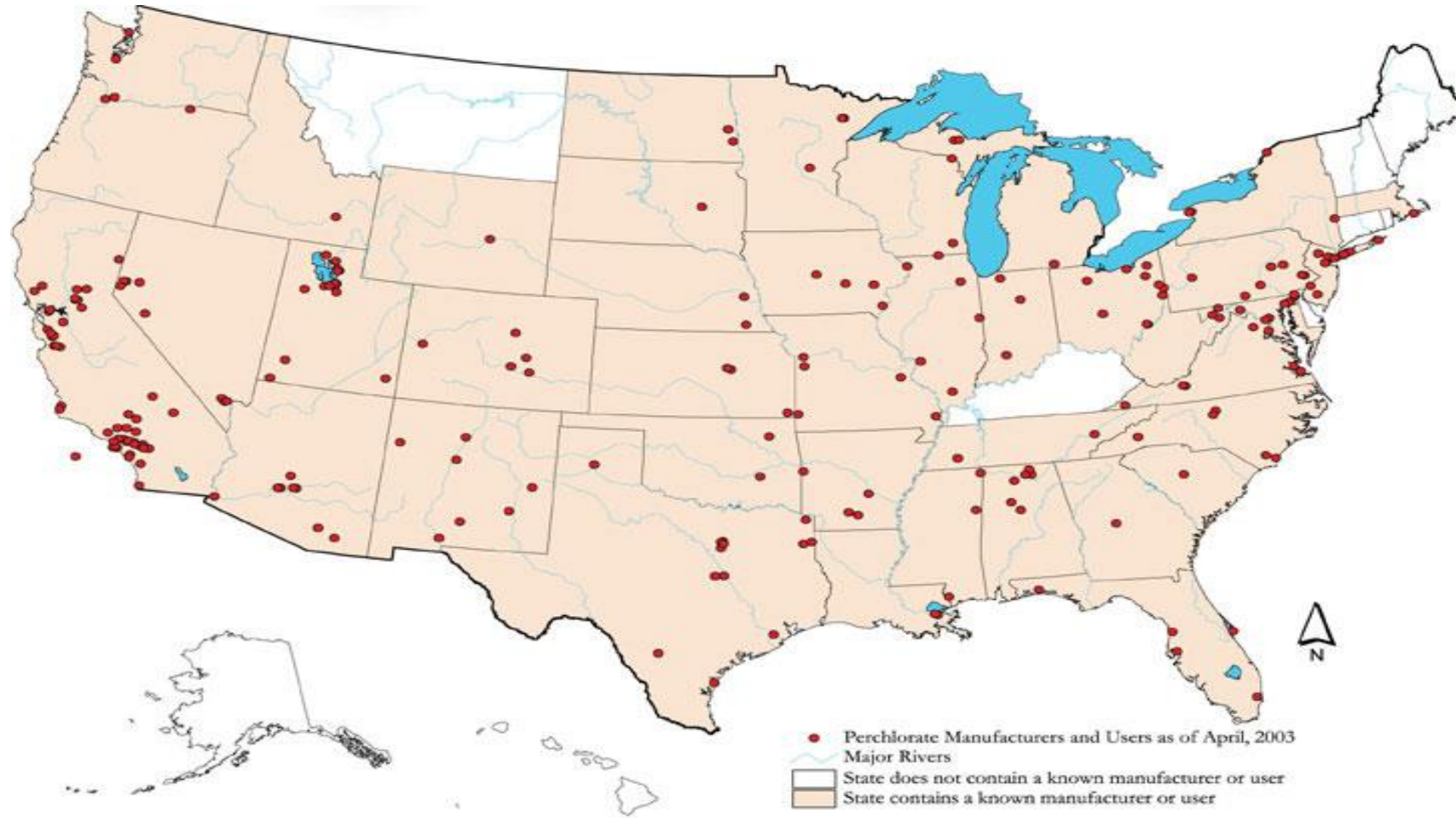

Map provided as public information on USEPA website - 2005

Possible Perchlorate Pathway

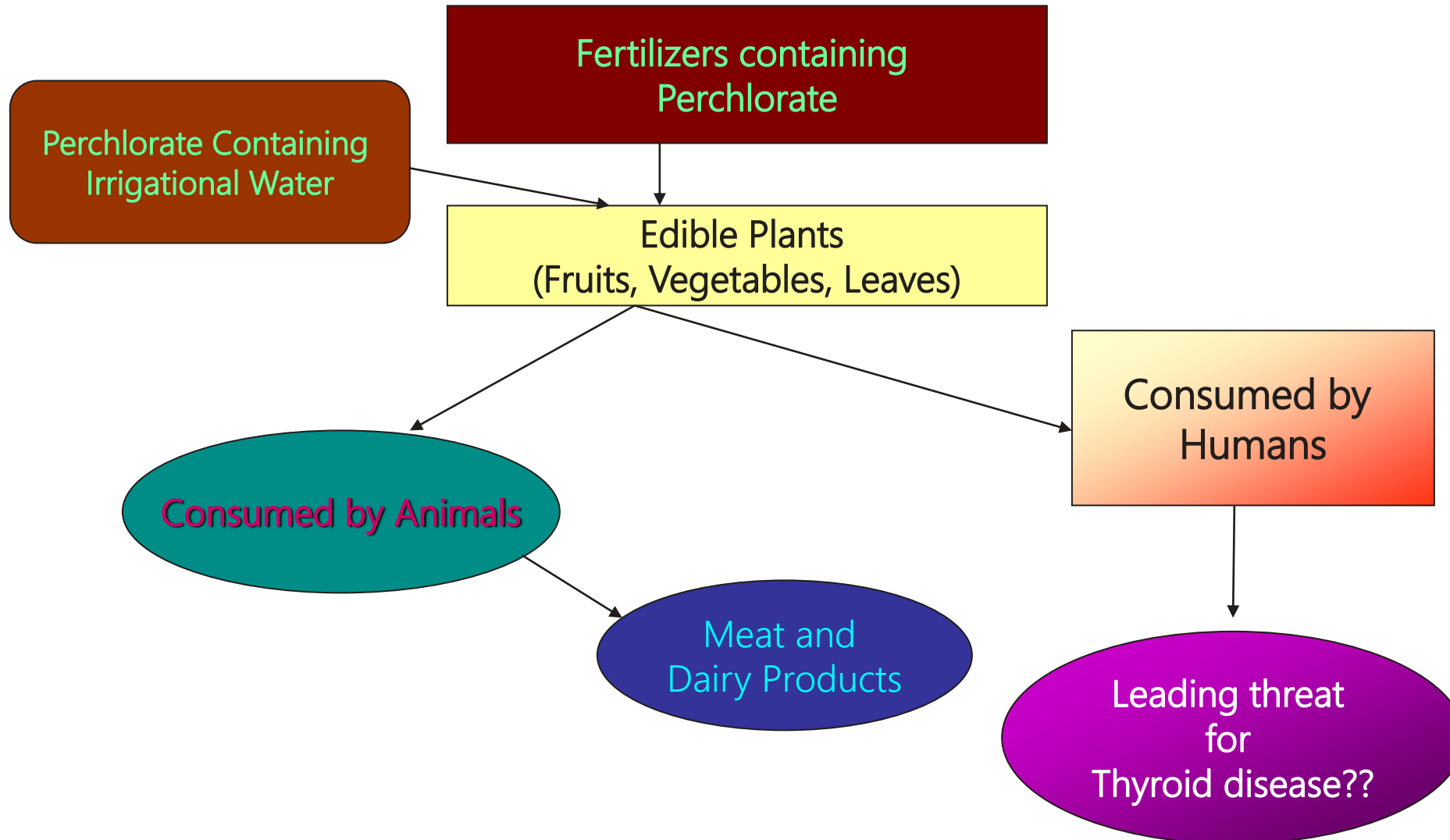

Regulatory Timelines

- Under SDWA, USEPA determines which chemical contaminant to regulate (CCL – Contaminant Candidate List)
- Perchlorate was part of
 - CCL1 – 1998
 - CCL2 – 2005
 - CCL3 – 2009
 - UCMR 1 – 2005
 - USEPA decision to regulate – Feb 2011
- In 2006 – State of Massachusetts regulates at 2ppb
- In 2007 – State of California promulgated at 6ppb
 - Most recent news – State of CA moving to regulate at 1ppb levels
- 12 Other states has adopted non-enforceable levels

Reference document: <https://www.regulations.gov/document?D=EPA-HQ-OW-2009-0297-0685>
<https://www.epa.gov/dwstandardsregulations/perchlorate-drinking-water>

USEPA Methods

•Perchlorate

- USEPA 314.0 enhanced
- USEPA 314.1
- USEPA 314.2
- USEPA 332.0 / SW846 6860
- USEPA 331.0 / SW846 6850

USEPA Methods

- Perchlorate

- USEPA 314.0 enhanced

- ~~•USEPA 314.1~~

- ~~•USEPA 314.2~~

- USEPA 332.0 / SW846 6860

- ~~•USEPA 331.0 / SW846 6850~~


US EPA method
314.0 enhanced
Perchlorate
by Suppressed Conductivity

USEPA method 314.0 (enhanced)

AW US6-0071

AW US6-0241


Using a Functionalized Monolithic Column or Anion Exchange Column

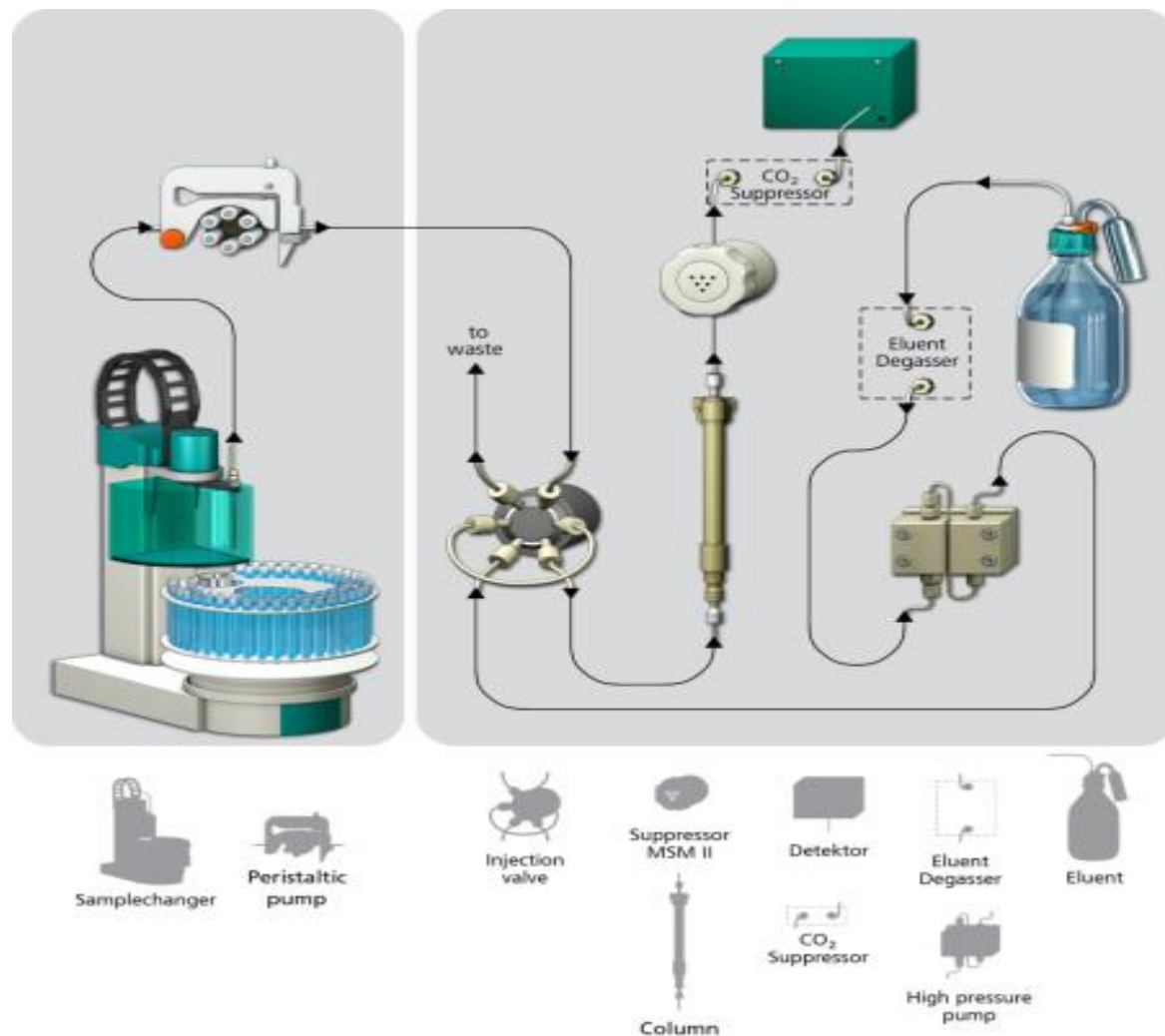

USEPA 314.0 - Perchlorate Standard in Matrix

Eluent flow	0.7 mL/min
Eluent	10.5mM Na ₂ CO ₃ + 25% Acetonitrile
Column	Metrosep ASupp7-250
Column temperature	45°C
Sample volume	1000 µL
Detector	Suppressed Conductivity

Component Name	Retention Time, mins	Concentration, mg/L
Chloride	7.5	1000
Sulfate	11.5	1000
Perchlorate	26.8	0.005

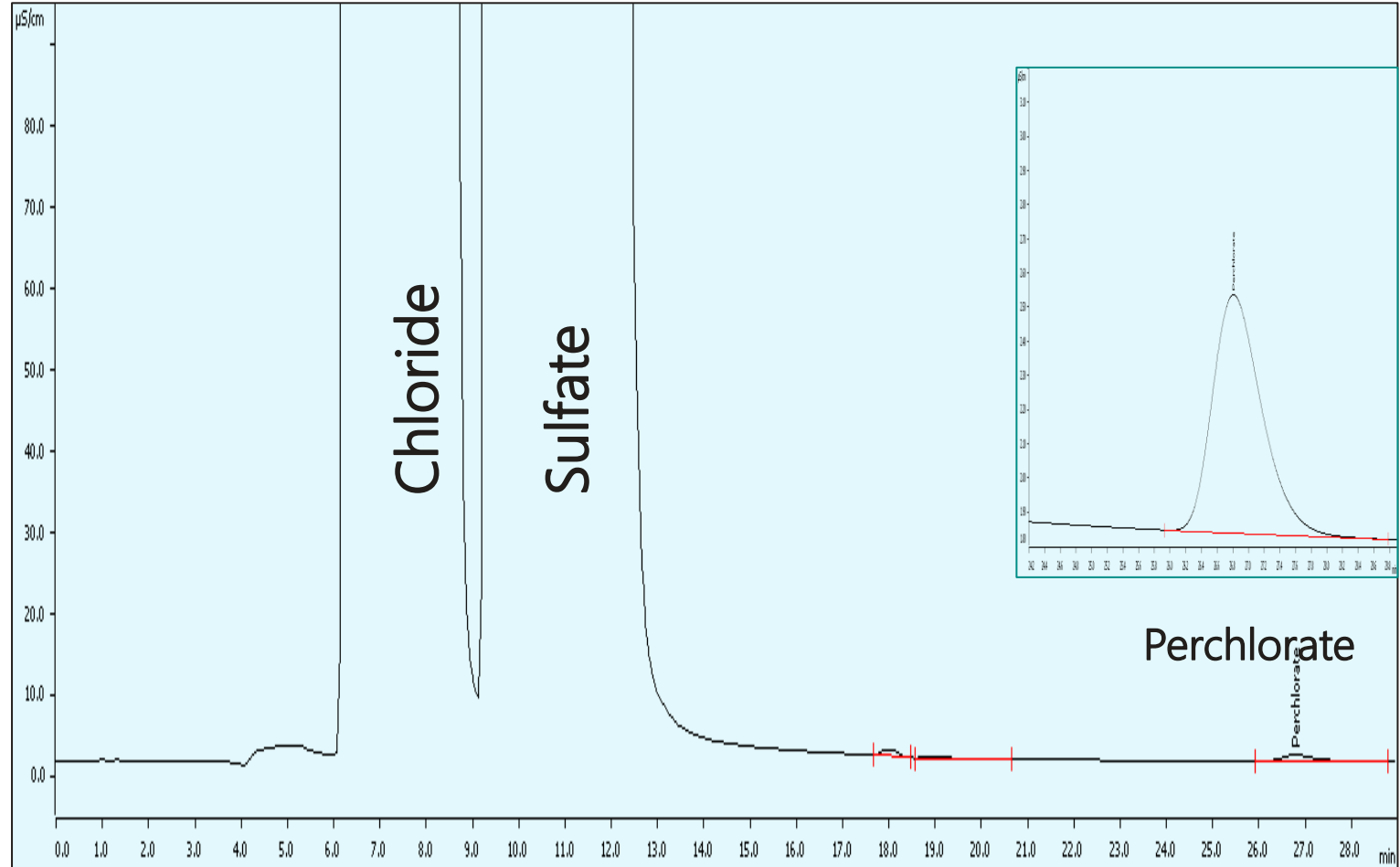


USEPA 314.0 MCT Study

Eluent flow	0.7 mL/min
Eluent	10.5mM Na ₂ CO ₃ + 25% Acetonitrile
Column	Metrosep ASupp7-250
Column temperature	45°C
Sample volume	1000 µL
Detector	Suppressed Conductivity

Component Name	Retention Time, mins	Concentration, mg/L
Chloride	9.28	50 - 1000
Sulfate	19.7	50 - 1000
Perchlorate	24.8	0.002


Maximum Conductivity Threshold Study (MCT Study)

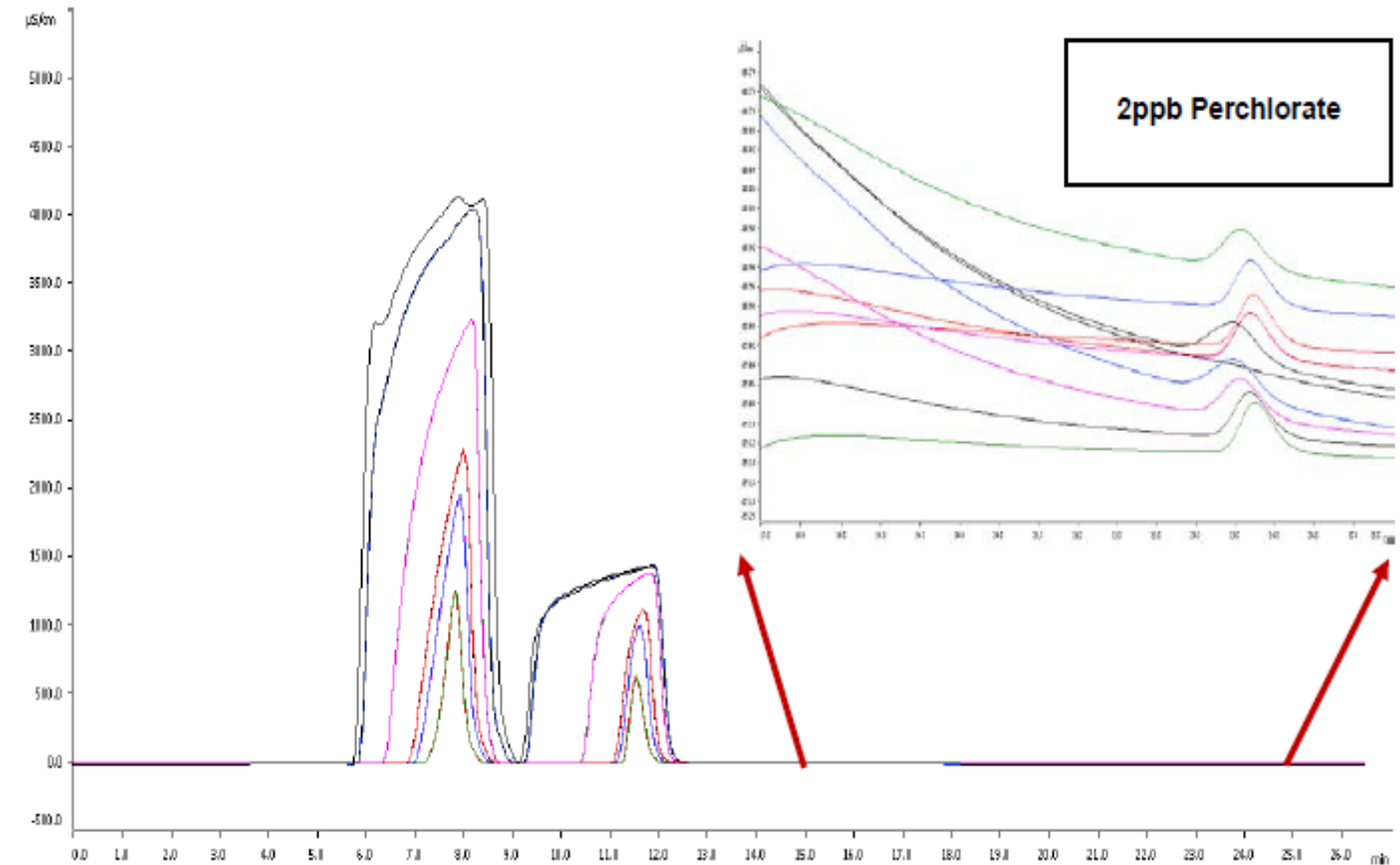

MCT Study Data (USEPA method 314.0, section 9)

MCT Study Data	
	Perchlorate, PGF
3000ppm TDS Matrix (n=5)	1.01
1500ppm TDS Matrix (n=5)	0.98
750ppm TDS Matrix (n=5)	0.94
300ppm TDS Matrix (n=5)	0.95
150ppm TDS Matrix (n=5)	0.93
50ppm TDS Matrix (n=5)	0.96

Major Interference – PCBSA

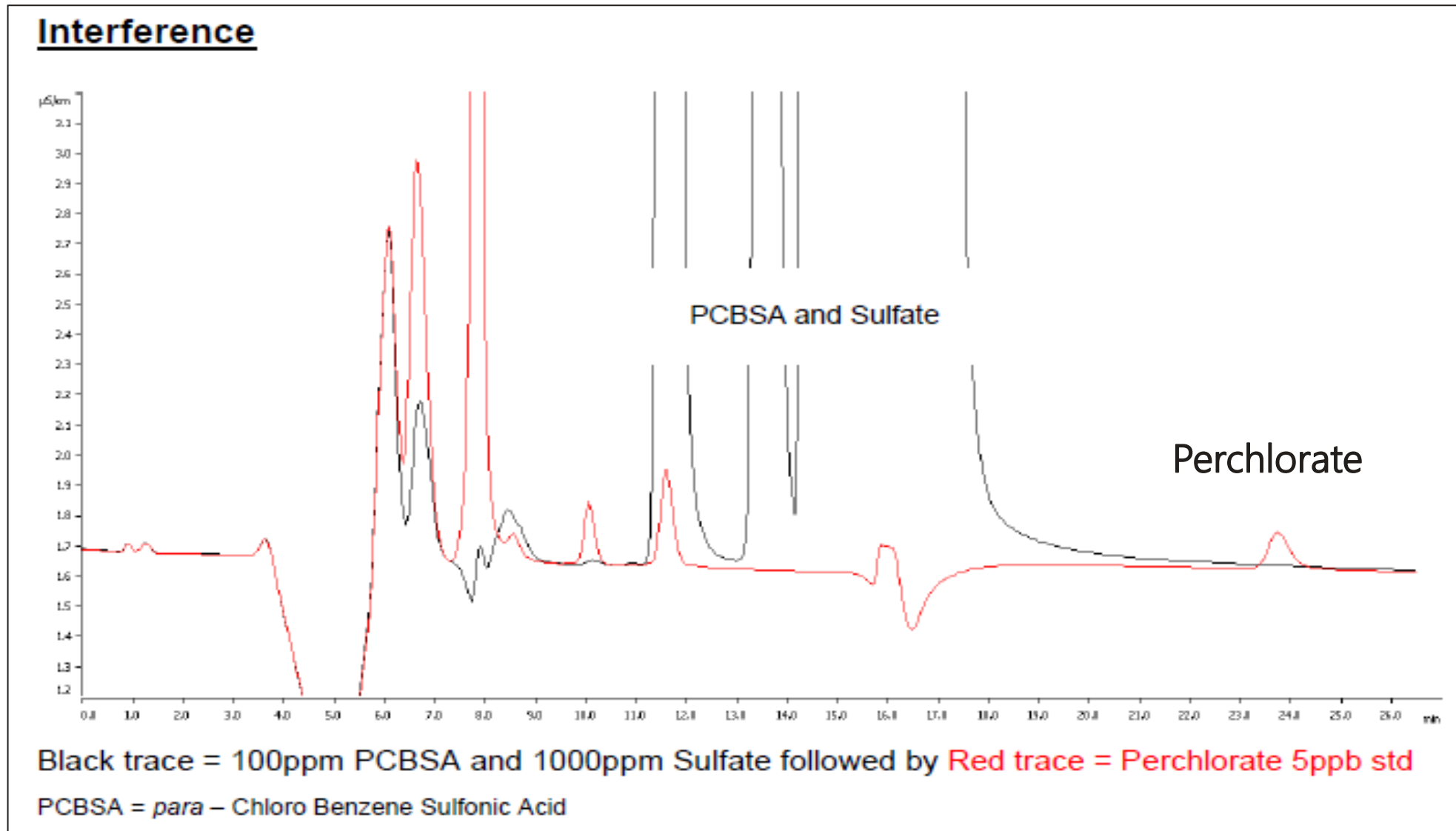

USEPA Method 332.0



+




OR




IC-MS

Electro Spray Ionization

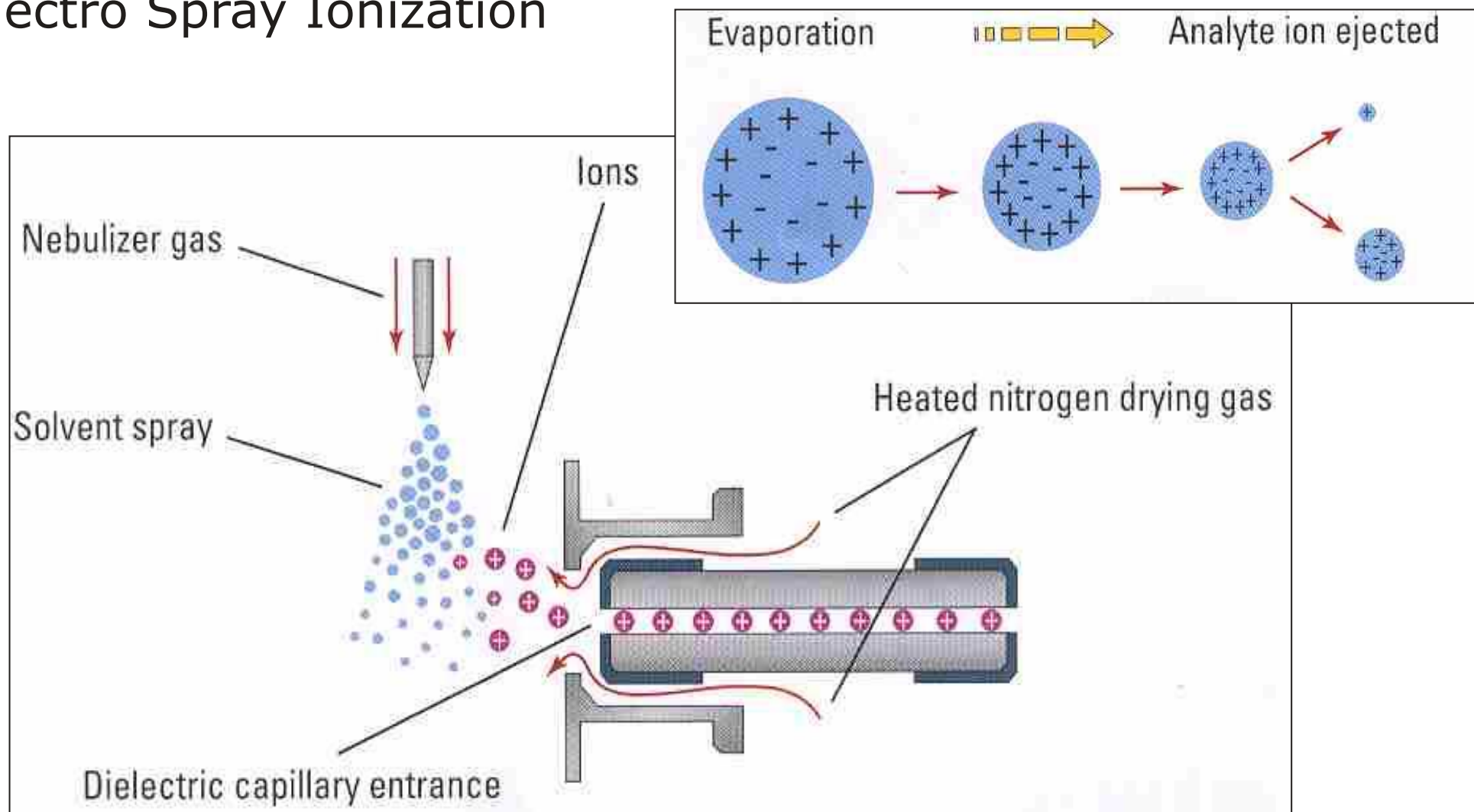

Perchlorate Mass

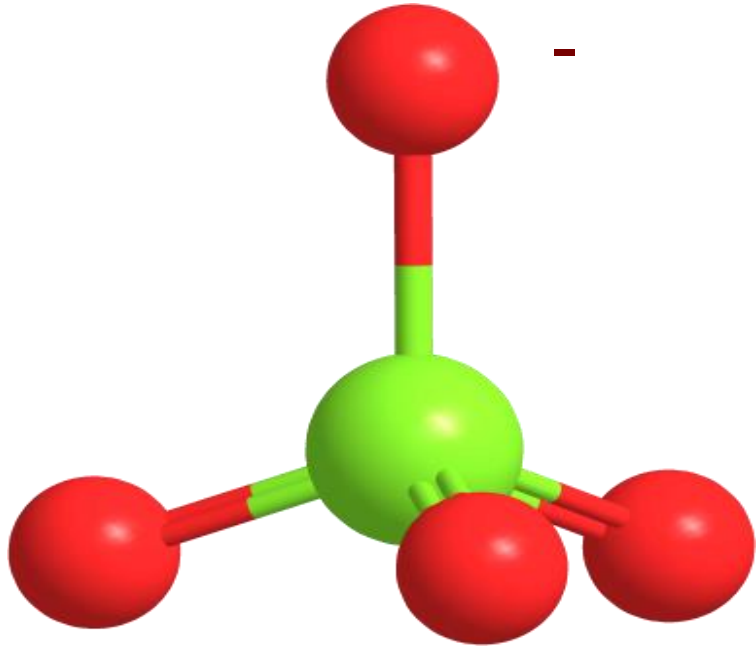

Chlorine (Cl) = 35

+

4 x Oxygen (O) = 4 x 16 = 64


Perchlorate mass (Cl³⁵) =
35 + 64 = 99

→ 75%

Perchlorate mass (Cl³⁷) =
37 + 64 = 101

→ 25%

Perchlorate Mass / interference

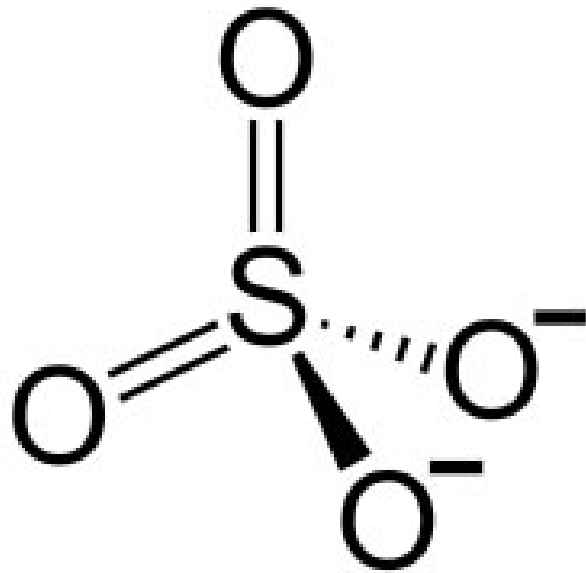

Sulfate Ion

Sulfur (S) = 32

+

4 x Oxygen (O) = 4 x 16 = 64

$$\text{HSO}_4^{-1} (\text{S}^{32}) = \\ 1 + 32 + 64 = 97$$

$$\text{HSO}_4^{-1} (\text{S}^{34}) = \\ 1 + 32 + 64 = 99$$

→ 4%

4% of 1000 parts per million = 40 parts per million OR
40,000 parts per billion

USEPA method
332.0 / SW846 6860
Perchlorate by ICMS / ICMSMS

In Collaboration with
USEPA (ODW-OGW / OSW)



Metrohm Advanced IC – Agilent iQMSD



Features

- Advanced Metrohm IC
 - Inline Cross flow filtration (optional)
 - Variable Volume Injection (optional)
- Agilent iQMSD
 - Standard ESI
- Open Lab CDS software for IC and MS control

Agilent Open Lab CDS



Instrument Status

Dashboard

858.0020 Professional IC Sample Processor

SQ

SQ

Idle

EMF ✓



858.0020 Professional IC Sample Pro...

Idle



Metrohm IC

Idle


Pressure MPa


Pressure MPa

Metrohm IC

Idle


55.0 °C




Metrohm IC

Idle


μS/cm


μS/cm

Metrohm IC

Idle


2


Last step
0.8 min

Metrohm IC

Idle



MS

A/S

IC

0.00 / 0.00


Instrument Idle



 On

 Off

USEPA method 332 / SW846 6860

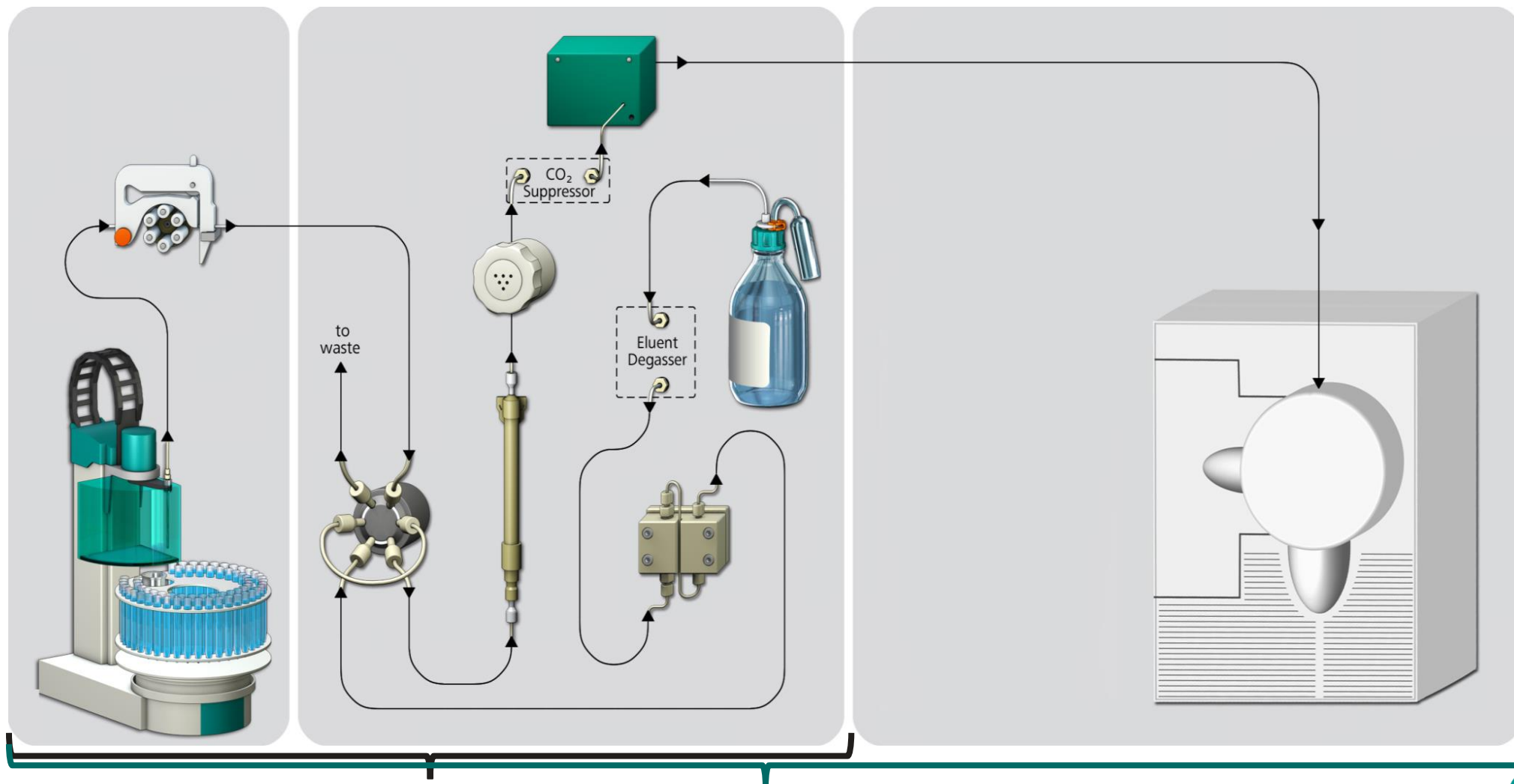

USEPA Method 314.0

USEPA Method 332.0

Chromatography

USEPA 332.0 –Calibration 0.5ppb & 25ppb - ICMS

Eluent flow	0.7 mL/min
Eluent	10.5mM Na ₂ CO ₃ + 25% Acetonitrile
Column	Metrosep ASupp7-250
Column temperature	45°C
Sample volume	100 µL
Detector	Agilent iQ MSD




Perchlorate




Perchlorate Calibration

AW IC US6-0240
AW IC US6-0244
AW IC US6-0291

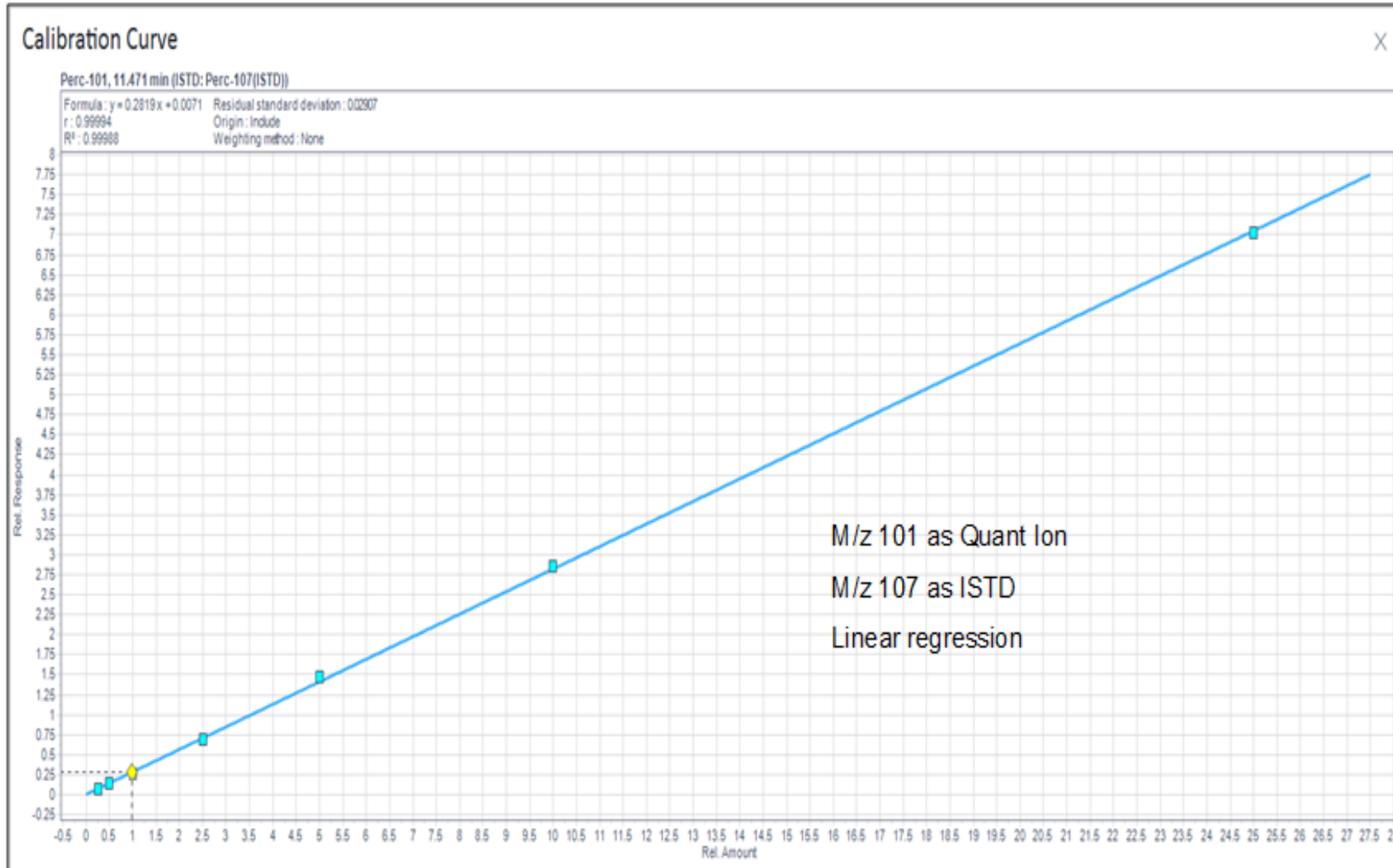


m/z 101
Range for
Calibration Standards
0.25ppb to 25ppb (ClO_4^{1-})

Real World Samples (example of Samples from CA)

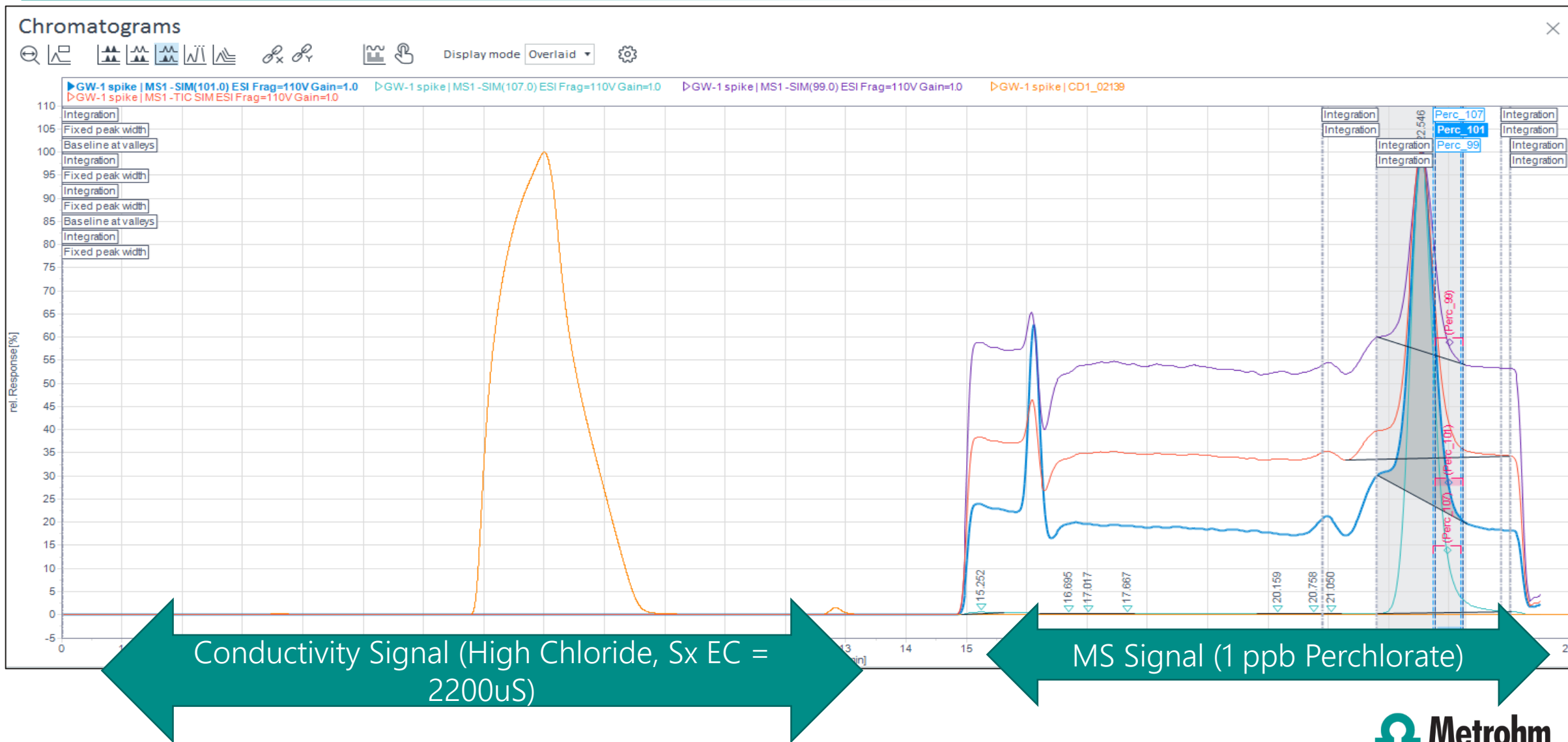

Real World Samples (example of Samples from Houston, TX)

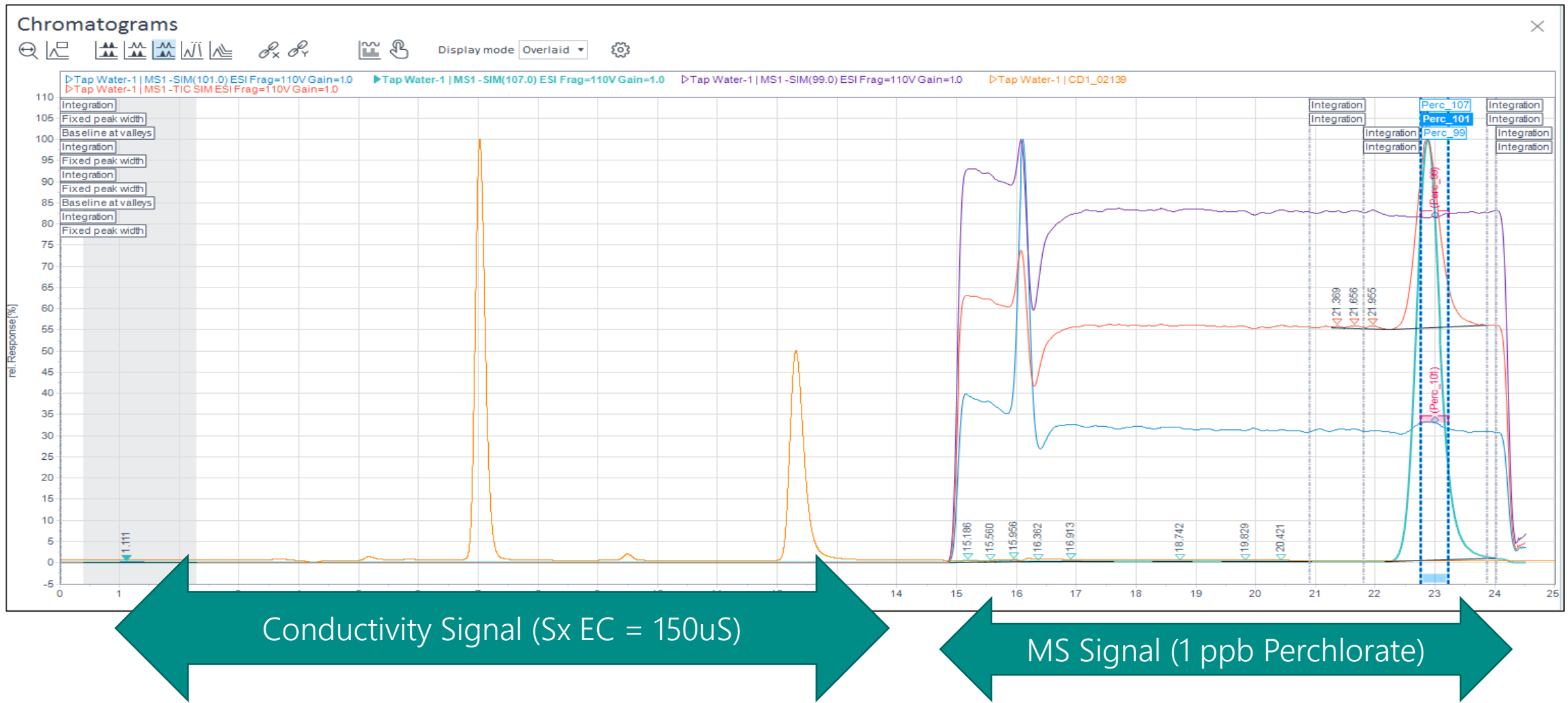

Reproducibility Data


Sample ID	m/z 99	m/z 101
0.5 µg/L standard	0.475	0.448
0.5 µg/L standard	0.468	0.472
0.5 µg/L standard	0.459	0.506
0.5 µg/L standard	0.463	0.483
0.5 µg/L standard	0.473	0.483
0.5 µg/L standard	0.475	0.495
0.5 µg/L standard	0.486	0.495
Average	0.4713	0.4831
Std. Dev	0.01784	0.01899
Calculated MDL (µg/L)	0.056	0.060

Sample ID	m/z 99	m/z 101
3000ppm Matrix + 1µg/L ClO ₄	0.96	1.08
3000ppm Matrix + 1µg/L ClO ₄	1.01	0.98
3000ppm Matrix + 1µg/L ClO ₄	1.01	0.99
3000ppm Matrix + 1µg/L ClO ₄	0.93	0.95
3000ppm Matrix + 1µg/L ClO ₄	1.08	0.96
3000ppm Matrix + 1µg/L ClO ₄	1.02	0.98
3000ppm Matrix + 1µg/L ClO ₄	0.97	0.92
Average	0.9971	0.9800
Std. Dev	0.04889	0.05000
Calculated MDL (µg/L)	0.154	0.157

What is Ion Suppression

- In LCMS / ICMS world, this is commonly known term
- Explanation of "ion suppression"


Effects of Ion Suppression

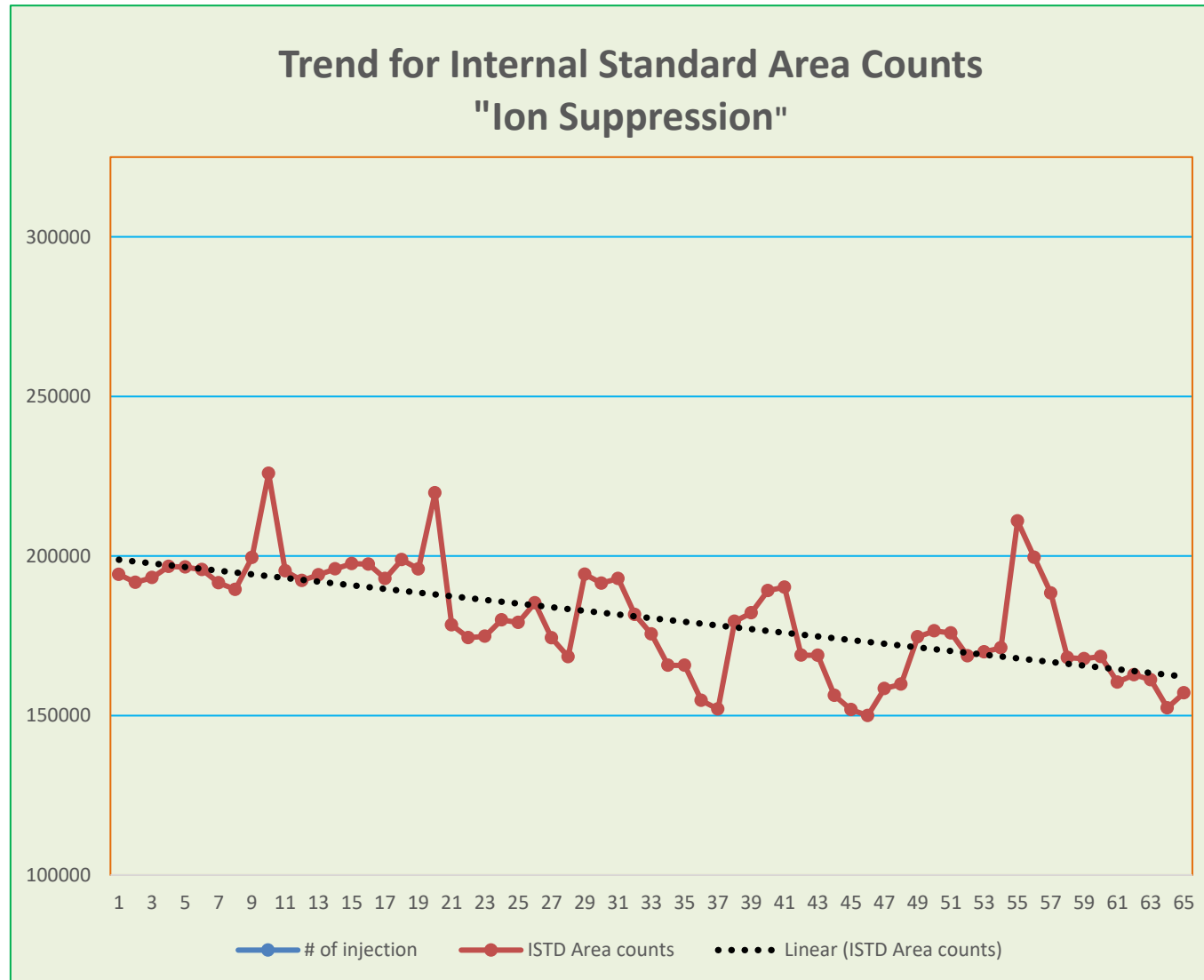

- Trendline indicates and demonstrates the "ionic suppression"
- However, "ionic suppression" has no significant loss in ESI performance



Before analysis

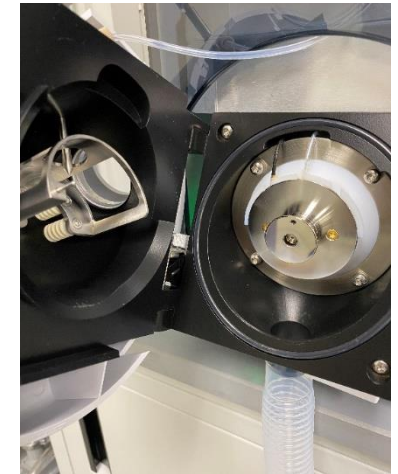

After 150+ injections

Summary

- Metrohm and Agilent Developed
 - Single software platform for ICMS analysis
 - Simple, rugged and isocratic analysis for Perchlorate
 - Environment friendly Carbonate chemistry for analytical column
 - Chromatographically resolved challenging matrices and interferences like
 - 3000 parts per million TDS (1000ppm each of Chloride, Carbonate and Sulfate)
 - Industrial surfactant like p-Chloro Benzene Sulfonic Acid (PCBSA)
- 100% Solvent compatible IC hardware allowed to hyphenate with
 - Single Quad MSD

Thank You

Be one of the forward thinking organizations that have experienced our full product portfolio