

Amperometric CN - Principles, Practice & Operational Benefits



A guide to selecting your CN method



Agenda



01 Amperometric CN - fundamentals

02 Amperometric CN - “practicals”

Performance

Method variants: Total CN, Available (WAD) CN, Free CN

Why use amperometric CN?

Instrumental considerations

03 Conclusions and questions

HELLO!



My name is Ilkka Lahdesmaki

Chief Scientist

You can find me at ilkka@flowinjection.com





1




AMPEROMETRIC CN FUNDAMENTALS

What is amperometry?

How does it measure cyanide?

Amperometry Principles

- Amperometry = measurement of electric current
- In chemistry, origin of current = electrochemical (redox) reaction
 - Requires **charge transfer**

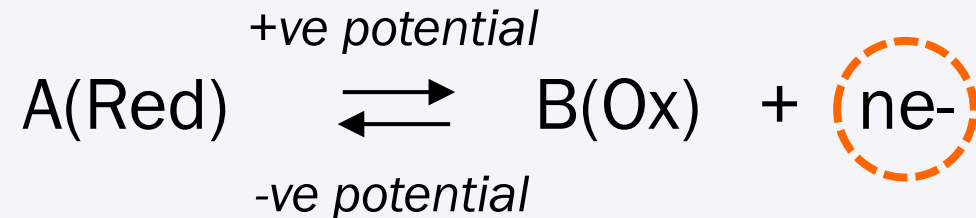


- OILRIG = Oxidation Is Loss, Reduction Is Gain (of electrons)
- Use a **potentiostat** to generate & measure current



Instrumentation & Methods


- Generation of amperometric signal relies on **mass transfer**
- Steady signal requires steadily moving solution
 - Can be implemented on a flow analyzer or using a stirred container
 - Cannot be run manually or on a discrete analyzer

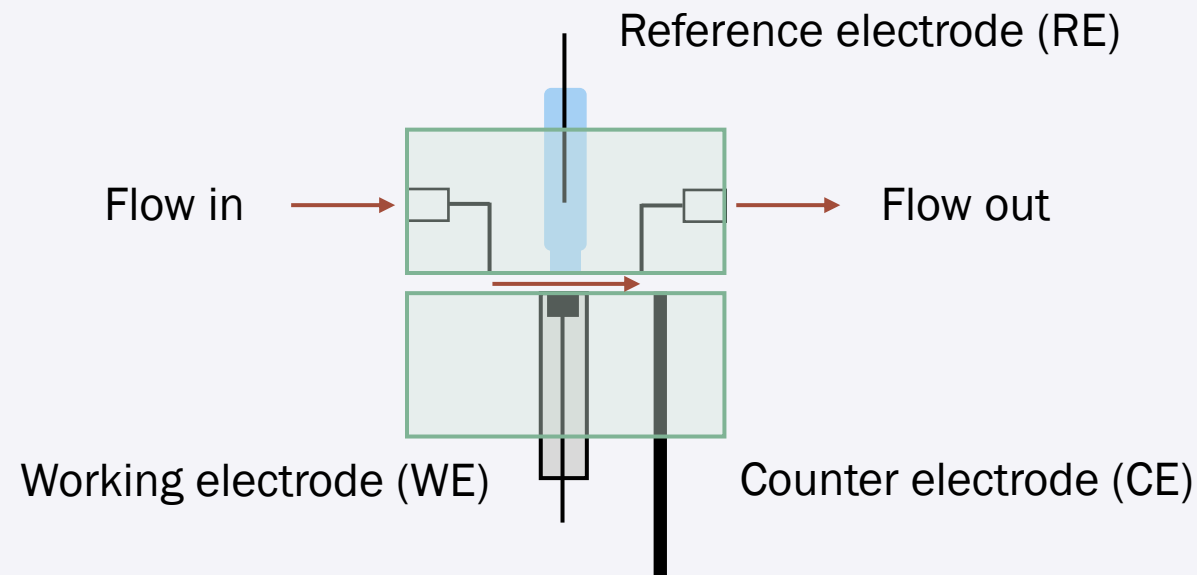

Instrumentation & Methods


◦ Potentiostat

- Applies a potential on working electrode (WE) to drive redox reaction
- Reference electrode (RE) is the reference point for that potential
- Measures current that flows through WE & counter electrode (CE)

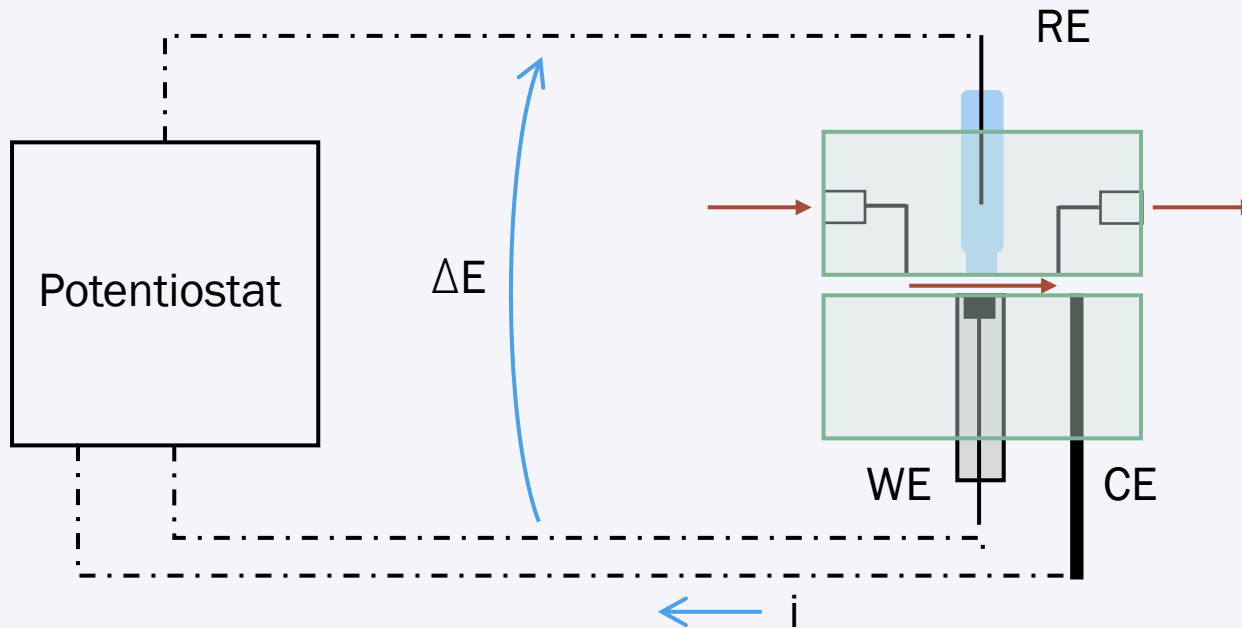


CN by Amperometry

- Detector configuration
 - WE = silver (Ag) electrode
 - RE = Ag/AgCl electrode
 - WE potential = 0.0 V

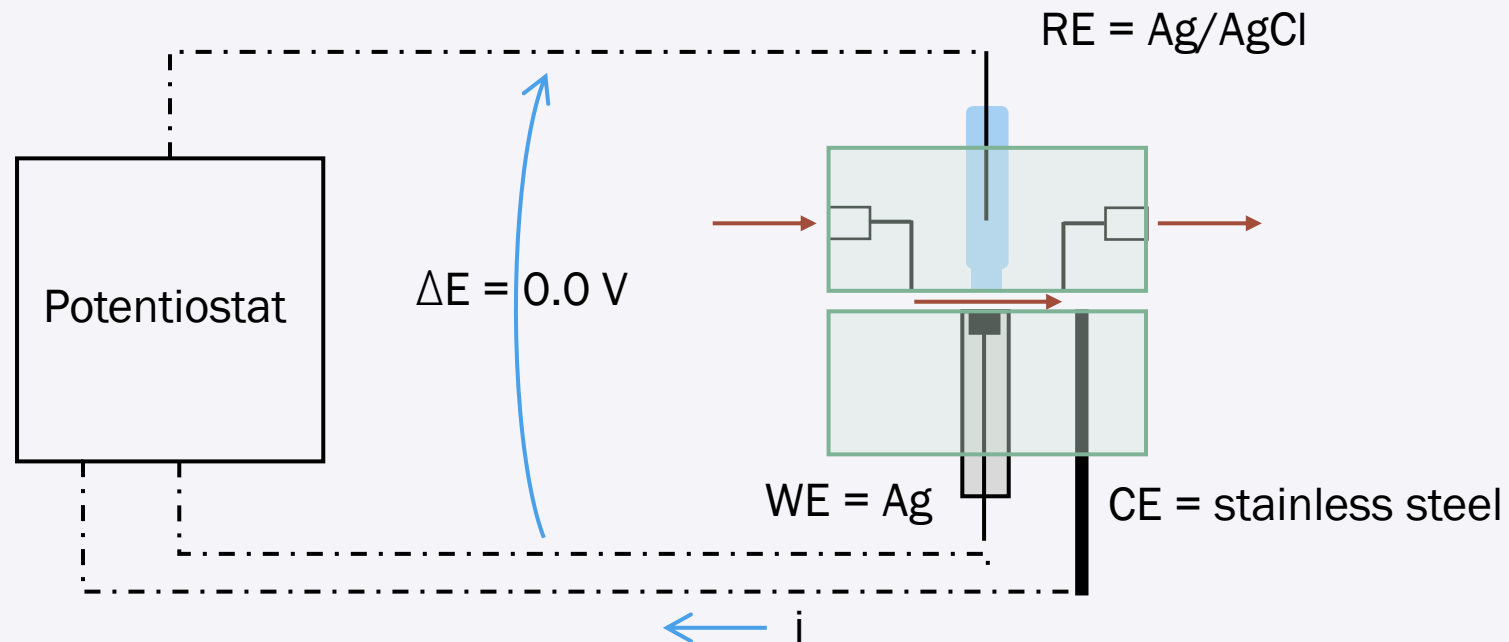


CN by Amperometry

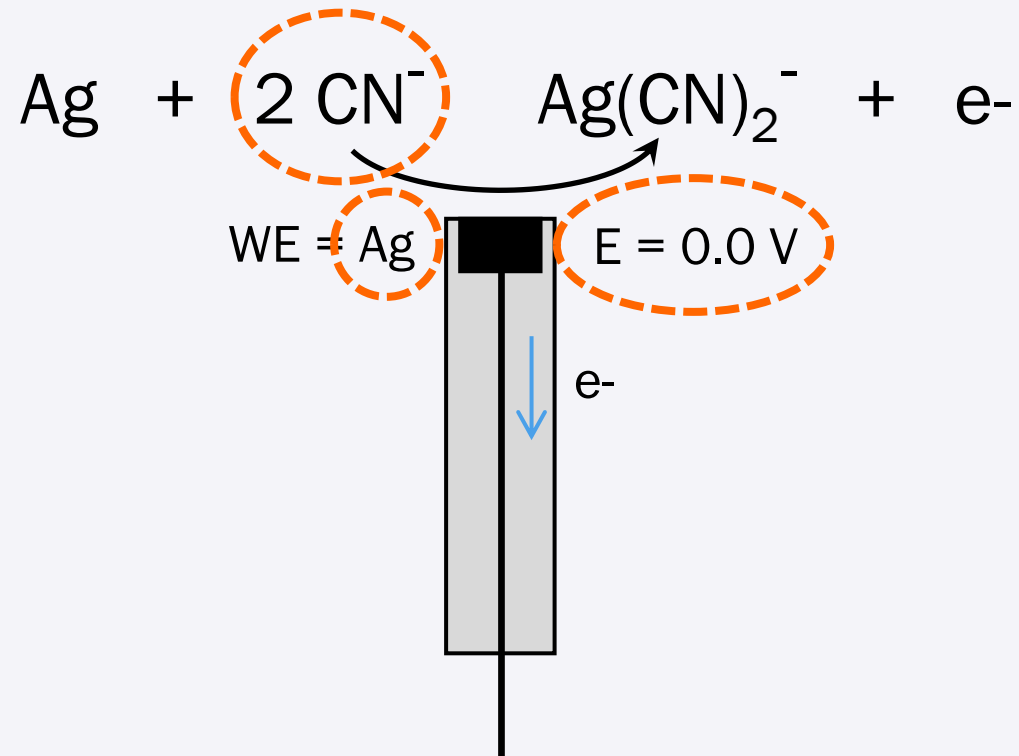

- CN⁻ reacts at Ag electrode, giving rise to a current (flow of e⁻)

CN by Amperometry

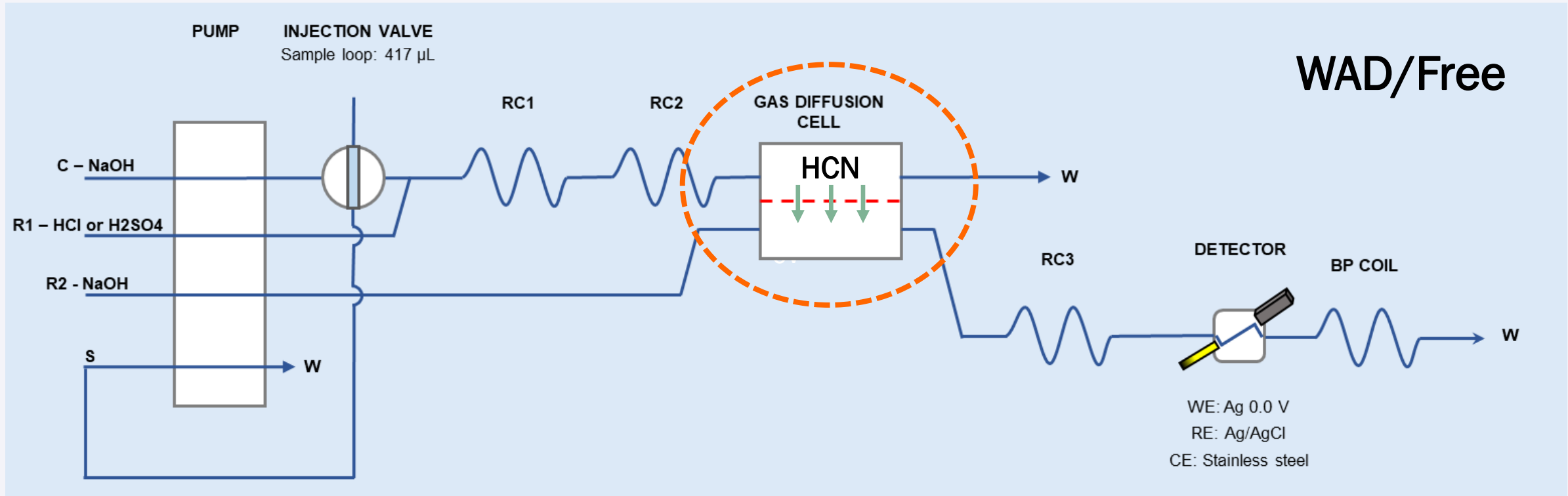

- All official amp CN methods rely on the concept of **gas diffusion**



AMPEROMETRIC CN “PRACTICALS”


How does amperometric CN perform?

Why use amperometric method for CN?


How to use amperometric method for CN?

Performance


- MDL: ~0.5 ppb (available/free), ~1 ppb (total)
- Range: usually up to 500 ppb, possible to adjust w/ sample loop size
- Throughput: 20-40 samples/h (per ASTM D7511-12)

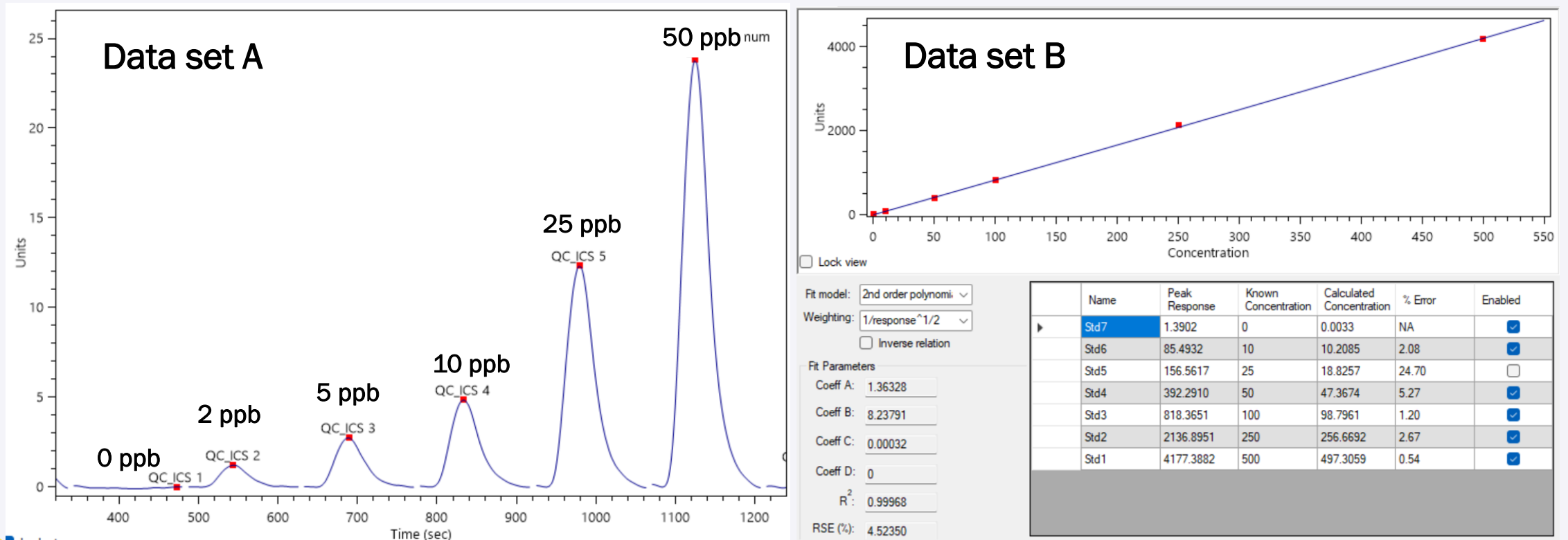



How to use amperometric CN?

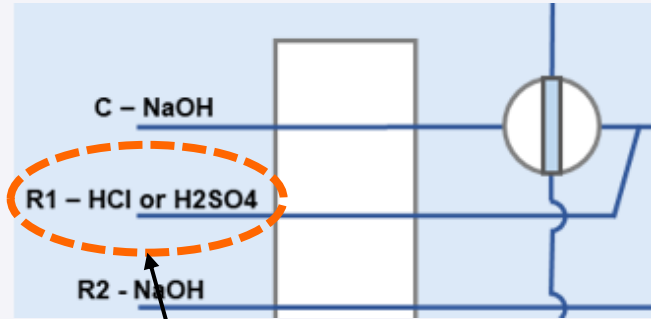


R1 formulation

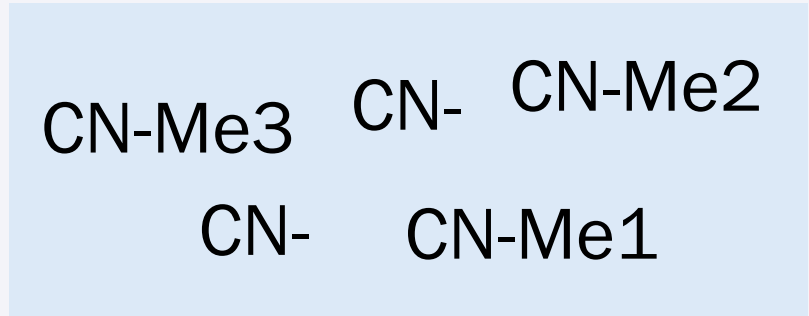


CN pool
in a sample

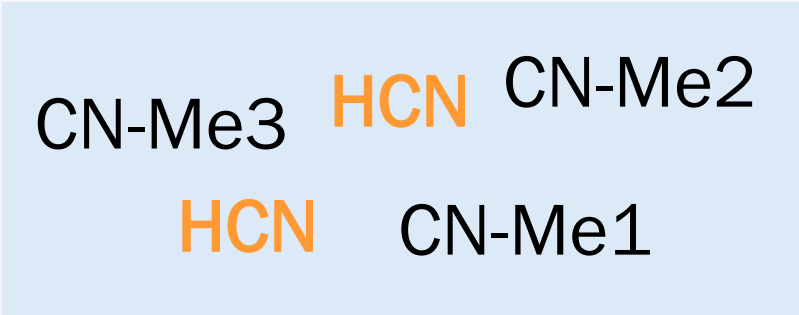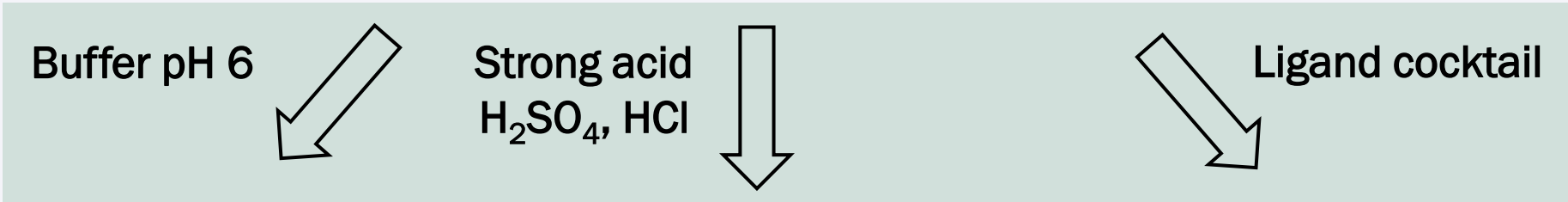

Free CN

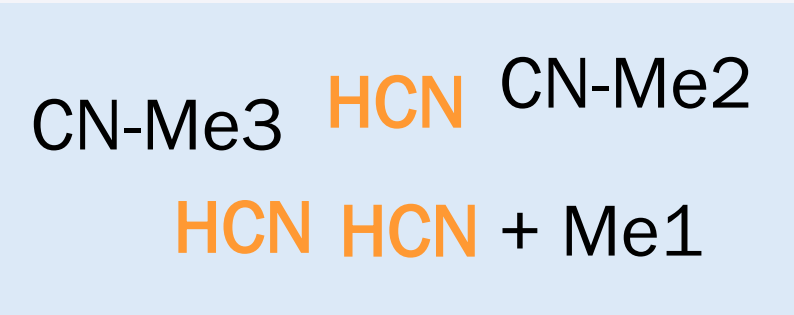

Available CN by acid dissociation
(a.k.a. WAD CN)

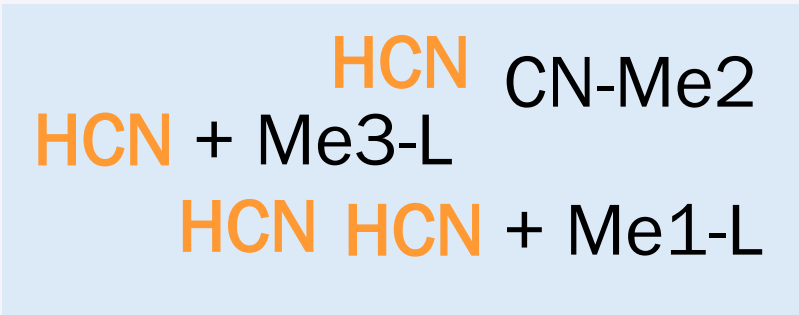


Available CN by ligand displacement
(a.k.a. WAD CN)



How to use amperometric CN?


CN pool
in a sample

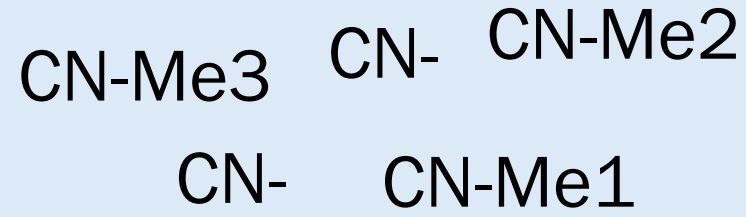

Strong acid
Reducing agent


↓

UV radiation

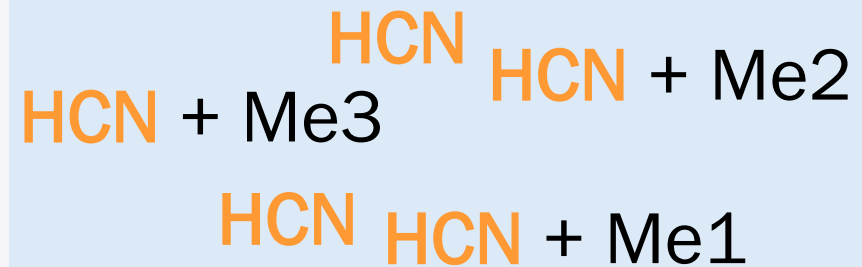


Total CN


CN by Amperometry

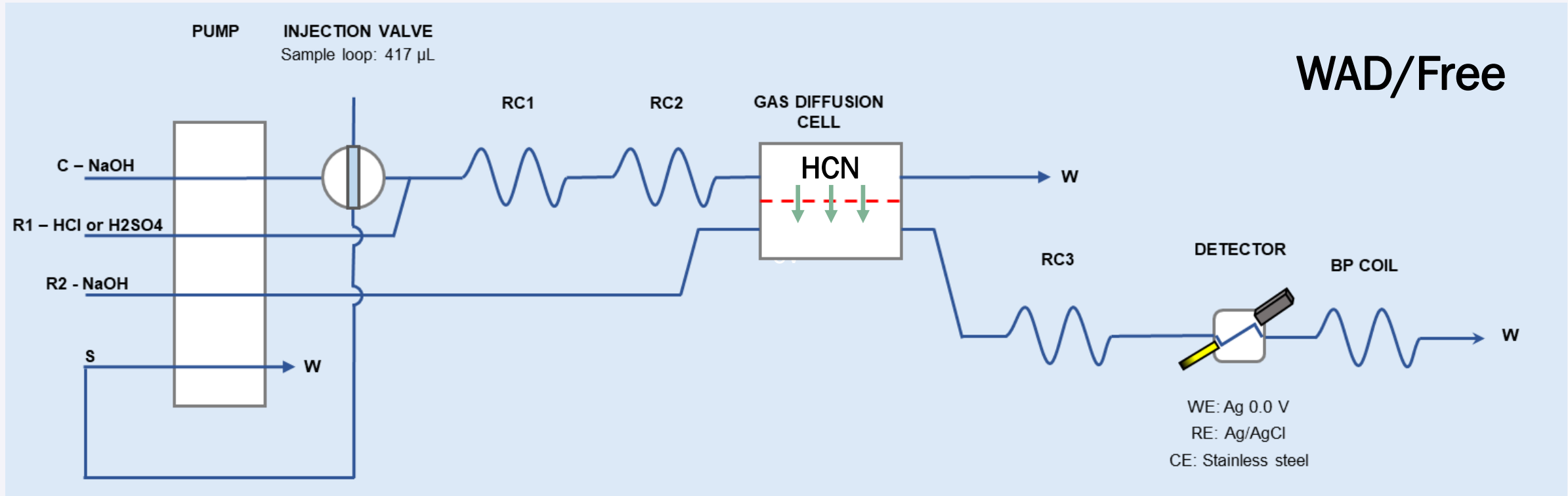

CN by Amperometry

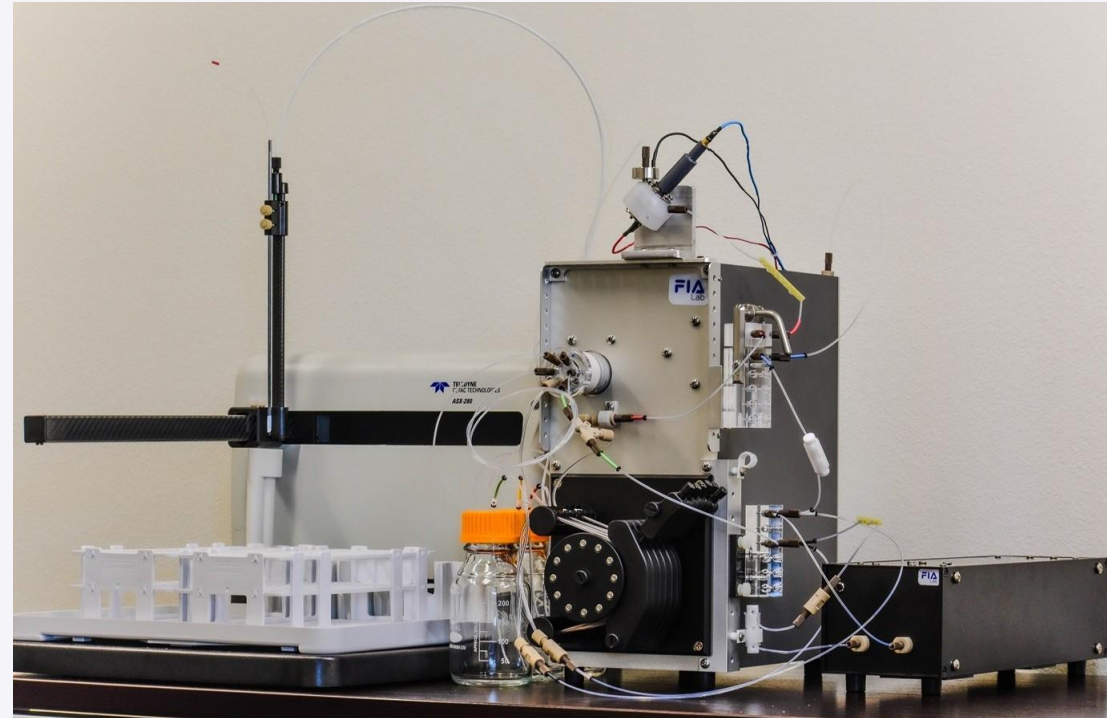

← ~10" →

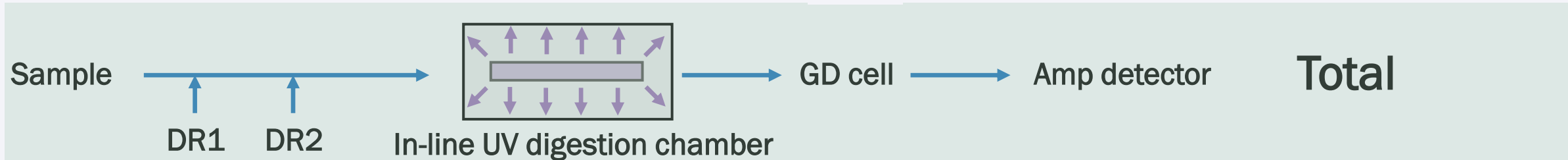

Why use amperometric CN?

- Simplicity
 - No need to deal with toxic chemicals (pyridine, Chloramine T)
 - Reagents simple to prepare, affordable (acid, base)
- Reliability
 - Distillation can result in false positives, UV digestion is less prone to that
- Approved methods available for different CN “classes”
 - Free CN
 - Available (a.k.a. WAD) CN (weak / intermediate CN-metal complexes)
 - Total CN
- Automated removal of sulfide interference



How to use amperometric CN?

- Many variant methods (* = included in 40 CFR Pt 136/141)

Free	• Free CN (buffer at pH 6): ASTM D7237* ¹³⁶ ; Uncomplexed CN
WAD	• Available CN (acid dissociable): EPA OIA-1677* ¹³⁶ ; Ag, Cd, Cu, Zn compl. • Available CN (acid dissociable): ASTM D6888* ¹³⁶ ; Ag, Cd, Cu, Zn compl.
WAD	• Available CN (ligand displacement): EPA OIA-1677* ^{136,141} ; + Hg, Ni • Available CN (ligand displacement): ASTM D6888* ^{136,141} ; + Hg, Ni
Total	• Total CN (following manual distillation): ASTM D7284* ¹³⁶ ; + Fe, Co, Au • Total CN (in-line digestion): ASTM D7511* ¹³⁶ ; + Fe, Co, Au

- “WAD” is not always clearly defined for available CN
 - Sometimes used for “acid dissociable”.
 - Sometimes used for “ligand displacement”.



Instrument considerations

- No two potentiostats are equal
 - Make sure the detector uses a model capable of determining ~1 ppb CN
- Find out whether the method implementations are practical
 - Avoid methods with “insane” sample segments, extreme low throughputs
- Instrument versatility
 - Easy conversion between Free, Available & Total CN setups
 - Get practical versatility with minimized capital investment
- Beware of heat in connection with UV digestion for Total CN
 - Heat can result in false positives (creates CN from SCN, CNO)
 - Can test by running ASTM D7511 “challenge matrix”
 - Some level of false positives invariable, “tolerance level” ~50 ppb





3



CONCLUSIONS

Conclusions

- Find out what data your customers need
 - Free, WAD, Total CN – or all of the above?
- What kind of sample load are you expecting for each?
 - Determines how many modules you need (1-channel / 2-channel)
- Talk to vendors of amp CN instrumentation
 - Practical MDLs, range
 - Ease of reconfiguration (1-channel used for WAD & Total)
 - Ease of operation
 - Customer support – do their technical staff understand amperometry?
 - If at all possible, talk to a customer reference





THANKS!



Any questions?

*You can find us at **Booth 51***

sales@flowinjection.com